

| | |
|---------|-------------------------------|
| 放大倍率: | 1x ~ 1000X. |
| 材質: | ABS 鏡片 PMMA |
| 電池容量: | 700mah. |
| 充電電池容量: | 2000mah. |
| 通訊功能: | WiFi Type-C Window Android |
| 照明: | LED. |
| 顯示屏: | 7 吋 IPS 觸控螢幕 |
| 顯示屏: | 8 吋 LED. |
| 顯示屏: | 可拆卸式顯示屏. |
| 認證: | CE FCC KC PSE |
| 重量: | 2.7 公斤. |
| 尺寸: | 25 * 25.5 * 21 公分 |





Explore the Micro World Around You



Bee wings



Ant



Bee leg



Longitudinal section
of corn seed



Slice ovary
of lily



Longitudinal section
of pumpkin stalks









How to Operate on



For Win10

Unpack and pick out microscope, connect the microscope and host computer via USB cable, search Camera in Windows for find it and click on.

(Notice: please see user manual for details)

For Win7 / Win8 / Mac

Step 1. Install software "HiView"

Step 2. Connect the microscope to PC via the USB cable, double-click to open the software

Step 3. Click "Device" in the menu bar, and then click "Open", select the microscope name "*****", and click "Open" to enter the preview screen

How to Use on Cellphone



Download the APP.



Long press the power button to turn on the microscope.



Turn on the phone's WIFI setting, find the name: "wifi_camera_wifi30_xxxx" and connect it.




Then you can click the APP and use



How to Use on Cellphone

 (Notice: please see user manual for details)

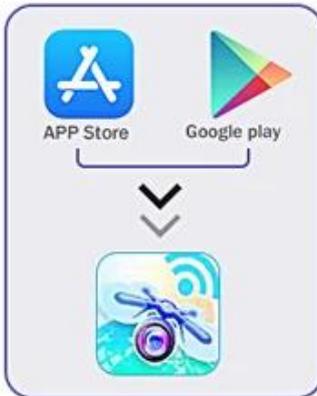


Step 1. Download the APP.



Step 2. Long press the power button to turn on the microscope.



Step 3. Turn on the phone's WIFI setting, find the name: "wifi_camera_wifi30_xxx" and connect it.

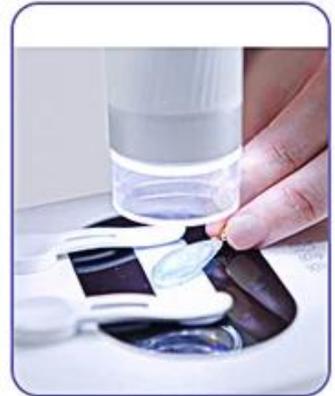

Step 4. Then you can click the APP

Wide Compatibility

Wi-Fi and USB 2 in 1 connection



iPhone



Android



Tablet



Mac




Windows



Wi-Fi



USB cable

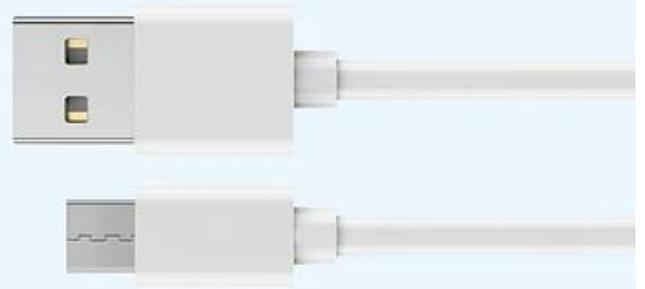



2700^{mAh}
High density
Li-ion battery

10^h
The base is in
continuous use

2^h
The monomer is
in continuous use



